

## Making Virtual, Reality

The GORE™ STA-PURE™ Bioprocess Configurator is an easy-to-use, interactive tool that enables users to work via the web to design and build customized fluid assemblies. This step-by-step online tool:

- Allows you to work independently to modify designs or interact directly with Gore engineers and authorized distributors
- Provides a flexible design mechanism via a drag-and-drop interface
- Reduces engineering design cycles and simplifies request for quote (RFQ) processes through an electronic approval system
- Enables 24-hours per day, 7-days per week design capabilities



# STA-PURE™ Bioprocess Configurator

*FOR CUSTOM FLUID ASSEMBLY DESIGN*



# STA-PURE<sup>SM</sup> Bioprocess Configurator

FOR CUSTOM FLUID ASSEMBLY DESIGN


## Customized Fluid Assembly Design Just Got Easier

The GORE<sup>SM</sup> STA-PURE<sup>SM</sup> Bioprocess Configurator enables the design of customized fluid transfer assemblies through our extensive portfolio of sampling devices, hoses and tubes, connectors, ends, peripherals and accessories.

Simply follow these steps to get started on a new configuration:

**One:** Go to [www.gore.com/pharmbio](http://www.gore.com/pharmbio), and click on the GORE<sup>SM</sup> STA-PURE<sup>SM</sup> Bioprocess Configurator link at the bottom, left-hand side of the screen. For first time users, click "Create an Account." Once you have created an account, you will be taken to the login page. Please login using your email address and password.

**Two:** To begin a design, simply drag and drop the desired product from the product palette onto the workspace. Save the drawing.

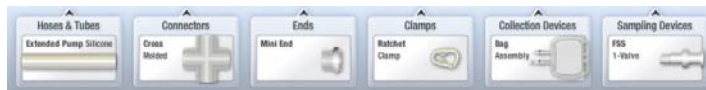

Once it is saved, you can modify the drawing before submitting it for a quote as well as during the request for quote (RFQ) process. Simply login to the system and submit changes. If you prefer, work with a Gore engineer or distributor partner to make any necessary changes.

**Three:** Submit the RFQ to Gore through the online system. An automated email will be sent to you and a Gore representative for review.

**Four:** Within three business days, a Gore representative will review the drawing and approve for quote or suggest alternative configurations. You will receive an email to approve the changes or make additional edits. Once approved, a Gore representative will email you a formal quote.

## Additional Features for Customizing the Design

**Properties:** The properties box automatically pops up when a new part is added. With this feature you can add measurements such as length, inner diameter, outer diameter, and backing hardware where applicable. The properties box provides a complete list of Gore products.



**Pressures:** Check this box to show the suggested maximum pressures for each part in the configuration.

**Metric or US Units:** For added convenience, design a configuration with metric or US units of measure.

**Custom Parts:** Check this box and custom parts within your design will be highlighted in yellow. Note: Custom parts may require additional lead times.

**Duplicate Selected Parts:** This feature reduces design time by enabling users to copy and paste parts that have already been designed. Click the part that you would like to duplicate and select this option, at the top right side of the screen.



## Contact Us

Visit [www.gore.com/pharmbio](http://www.gore.com/pharmbio) for more information on the GORE<sup>SM</sup> STA-PURE<sup>SM</sup> Bioprocess Configurator. For additional questions, contact a Gore representative or an authorized Gore distributor. If you are interested in receiving additional training on this tool, please contact a Gore representative.

To the best of our knowledge, the information contained herein is accurate. Final determination of the suitability of any information or material for the use contemplated and the manner of use is the sole responsibility of the user. The above information gives typical properties only.

GORE and design are trademarks of W. L. Gore & Associates, Inc. STA-PURE is a service mark of W. L. Gore & Associates, Inc. All rights reserved. © 2008 W. L. Gore & Associates, Inc. Printed in USA.

### W. L. GORE & ASSOCIATES, INC. | GORE PHARMBIO PRODUCTS

402 Vieve's Way • Elkton, MD 21921 • USA

Phone: 410.506.7787 • Toll-Free: 800.294.4673 • [pharmbio@wlgore.com](mailto:pharmbio@wlgore.com)

[gore.com/pharmbio](http://gore.com/pharmbio)

### DISTRIBUTOR INFORMATION

